

# CAGEWASH 2

## NEUTRAL DETERGENT CLEANER

### A NEUTRAL LOW-FOAM CLEANER FOR WASHER DISINFECTORS & MANUAL APPLICATIONS

PH neutral detergent which is ideal for use within cage washers, bottle washers and rack washers for removing organic soiling on all types of caging and associated equipment. It is also effective at inhibiting surface scale and mineral deposit build up.

Cage Wash 2 is economical, versatile and has multi-surface use, suitable on both hard and soft water applications. It prevents & removes scale build-up and is ideally suited to hard water due to its scale removing properties and excellent rinsing capabilities.

Suitable for use on low temperature washing below 50°C.

## Application

1. Place feed tube into bottom of container and check automatic feed equipment is switched on.
2. Inject at 4ml per 1 litre of water at a working temperature of 45 – 60 °C.

## Material Compatibility

Cagewash 2 is compatible with the following materials: Stainless Steel, Aluminium, Anodised aluminium, Chromium plated material, Synthetic materials, Rubber, Latex, Glass, Ceramics, Copper, Brass, Metal alloys.

## Warnings and Cautions

- Do not add any other chemicals, such as bleach or detergents. Added chemicals may destroy the effectiveness of this product.
- Wear goggles, plastic apron, rubber gloves, when changing containers.
- Consult the manufacturers recommended reprocessing instructions to determine cleaning agent compatibility and any additional procedures necessary to clean their particular medical device.
- Store in a cool place between 5 and 25°C.
- Dispose in accordance with local authority guidelines.



Available in 5L, 10L, 20L and 200L.