

# CAGEWASH 3

## ALKALINE DETERGENT CLEANER

### CONCENTRATED ALKALINE DETERGENT CLEANER

Heavy duty alkali detergent which is ideal for use on cage washers, bottle washers and rack washers for removal of organic & greasy soiling from caging and associated equipment. Safe to use on glass, stainless steel and plastics.

Cage Wash 3 is economical, versatile and has multi-surface use, suitable on both hard and soft water applications. It effectively removes heavy, baked on organic and greasy deposits often seen on larger caging and racks and has excellent rinsing capabilities.

Suitable for use on low temperature washing below 50°C.

## Application

1. Place feed tube into bottom of container and check automatic feed equipment is switched on.
2. Inject at 4ml per 1 litre of water at a working temperature of 45 – 60 °C.

## Material Compatibility

Cagewash 3 is compatible with the following materials: Stainless Steel, Aluminium, Anodised aluminium, Chromium plated material, Synthetic materials, Rubber, Latex, Glass, Ceramics, Copper, Brass, Metal alloys.

## Warnings and Cautions

- Do not add any other chemicals, such as bleach or detergents. Added chemicals may destroy the effectiveness of this product.
- Wear goggles, plastic apron, rubber gloves, when changing containers.
- Consult the manufacturers recommended reprocessing instructions to determine cleaning agent compatibility and any additional procedures necessary to clean their particular medical device.
- Store in a cool place between 5 and 25°C.
- Dispose in accordance with local authority guidelines.



Available in 5L, 10L, 20L and 200L.